**LJ CREATE**  
Learning for life

# Datalog Software

Quick Start Guide  
Guía rápida

## System Requirements

<b>Operating Systems:</b>	Windows XP or Vista
<b>Display:</b>	Minimum 800 × 600 with 256 colors Recommended: 1024 × 768 or above
<b>Minimum Software:</b>	Microsoft Internet Explorer 6.0 Windows Installer 3.1* Microsoft .NET Framework 2*

\* If Microsoft .NET Framework 2 or Installer 3.1 is not detected during the installation set up process, they will be automatically installed prior to the dataLOG application install. This will extend the installation time significantly.

### The dataLOG Software CD-ROM

A Windows installer package (**.msi**) of the dataLOG software is supplied on the CD-ROM.

The Drivers folder on the CD contains the USB device drivers for all models of the dataLOG loggers and the AMC232USBDH USB Serial Adapter cable. The dataLOG drivers (XP only) will install to the local machine at the end of the dataLOG software install.

A PDF format of the dataLOG Software Manual and a copy of this Quick Start Guide, are available in the PDF folder on the dataLOG CD-ROM. It is advised that the latest version of Adobe Acrobat reader software is installed to read these documents.

## Step 1: Install the dataLOG software

**Windows XP and Vista users:** Log on as ADMINISTRATOR (or with equivalent access rights) on the system.

**USB users:** Do **not** connect the dataLOG logger to the USB port prior to, or while, the dataLOG software is being installed.

**Current dataLOG users:** If you have an earlier version of dataLOG software on your computer, remove it before commencing the installation, i.e. use Add or Remove programs accessed via Start ▶ Settings ▶ Control Panel. You may need to restart your machine for configuration changes to be made.

**All users:** Insert the dataLOG CD-ROM into the CD drive. This is an auto play CD-ROM, meaning that Windows should detect the presence of the CD and automatically start up the dataLOG auto run program. If the auto run program fails to start, use Windows Explorer to navigate the CD-ROM contents and double click on the file **setup.exe**.

**Note:** *If .NET Framework 2 is not detected, it will now install – our experience is that this may take between 3 – 15 minutes.*

- Follow the on-screen instructions to complete the installation of dataLOG and its associated files onto the hard disk.

**Windows XP users:** The dataLOG (EasySense) device drivers will be pre-installed to the local drive at the end of the install. A message may appear for each device driver stating that they have not passed Windows Logo signing, please select 'Continue Anyway'. This message may appear more than once. When the installation is complete, eject the CD-ROM and store it safely.

- Restart your system.

## Step 2: Connect an Interface or Data logger to the computer

### Serial connection:

- Some data loggers, for example dataLOG 100, can connect to the computer via the serial port on the computer. If you use a serial port connected data logger note the serial port you have connected to e.g. COM 1, as this information will be required when you start the dataLOG software.

### USB connection:

**Windows XP and Vista users:** Log on as Administrator (or with equivalent access rights) on the system.

The first time a dataLOG unit is connected to a USB port on the computer, the USB drivers will need to be installed.

- Ensure the data logger is awake by pressing any of the buttons on the data logger. Connect the dataLOG unit to the USB port on the computer. Windows will automatically detect that a new device has been connected and a Wizard will be generated to aid installation.

**Window Vista users:** Select 'Locate and install driver software (recommended)'. You may be asked for permission to continue. If the PC is connected to the Internet, Windows Vista may try to find the driver automatically from Microsoft's Windows Update servers. At this point it will fail to find the drivers and you will be asked to install the drivers from the CD-ROM, insert the CD-ROM into the drive and click Next. You may be warned that the drivers have not passed Windows Logo signing, select 'Install the driver software anyway' at this point. The USB drivers are also copied to the local drive [default location is C:\Program Files\DHG\Drivers\USB Drivers folder] and can be installed from here if the CD-ROM can not be located.

**Windows XP users:** If prompted to connect to the Internet to find the drivers automatically from Microsoft's Windows Update servers select 'No, not at this time' and click on Next. Select to 'Install the software automatically'. You may be warned that the drivers have not passed Windows Logo signing, please 'continue anyway' at this point.

- Connect the dataLOG unit to every USB port on the computer so that the drivers can be installed for each port.

## Step 3: Start the dataLOG software for the first time

Start the dataLOG software e.g. Start ► Programs ► LJ Create ► dataLOG and click on dataLOG software.

## Step 4: Select a User level

The first time the dataLOG program is opened, a 'Select Program Level window will automatically open. Select a suitable user level.


- Level 1 is aimed at students up to 9 years old
- Level 2 is aimed at students between 9 – 15 years old
- Level 3 is aimed at students over 15 years old

**Note:** The program level can easily be altered at anytime using the Level icon once a logging option has been opened. Click on the Level icon and cycle through the levels. The currently selected level will be saved when the dataLOG software is closed.

## Step 5: Select an Interface and Connection

The first time the dataLOG software is opened, a 'Select Interface' window will automatically open to allow the correct interface and connection to be selected.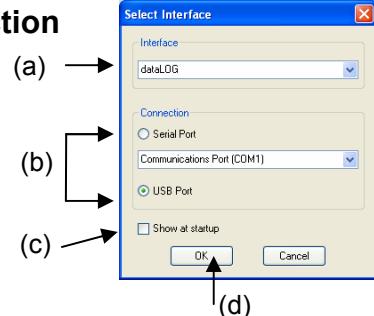

## English

- (a) Select the correct interface from the drop-down menu.
- **dataLOG** for a dataLOG type data logger.
  - **Example Data** for stored example data to show how sensor readings are displayed. If this option is selected you will not be able to collect data from your data logger.
- (b) Select the appropriate connection i.e. either Serial or USB Port.
- If you are connecting via a **USB** lead, select USB Port.
  - If you are using a **serial** lead, select Serial Port and then the number of the communications port used from the drop-down menu e.g. Communications Port (COM2).
  - If you are using a **USB/serial adapter**, select Serial Port and then the adapter from the serial ports drop-down list, for example, AMC232USBDH USB Serial Port (COM4).
  - If you are using **Bluetooth**, select Serial Port and then the serial port configured for Bluetooth from the drop down list, for example, Bluetooth Communications Port (COM4).
- (c) If these settings will stay the same, leave the ‘Show at startup’ box unchecked. If the box is checked (enabled) the ‘Select Interface’ window will automatically open when you next start the dataLOG software, so you can see if the currently selected Interface and method of connection are correct. This is useful if you use different methods of connection (for example, using both the dataLOG 100 type unit which connects via a serial port and dataLOG 120 type unit which connects via a USB port).
- (d) Click on OK. The program will save the current selection.

**Note:** To change this selection at any time, CANCEL any ‘failure to connect’ errors, select **Interface** from the **File** menu and alter the interface selection.

Once a connection is established the Home page will open, click on the appropriate icon.

## The Home page


Home is the main navigation page, which contains the icons for all available options.

The appearance of the Home page and the complexity of the toolbars and functions in the experiment options are altered by the User level selected. The more bars, the higher the user level.

The Home page at Level 1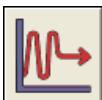

### EasyLog

The Graph window will open pre-set to record sensor values **continuously** as a line graph until stopped.


Click on the **Start/Stop** icon  to start recording. Logging will start with a time span of 30 seconds and when it has elapsed the time span will double automatically. This will continue until the recording is stopped by clicking on the **Start/Stop** icon. 

### Meters

The Meters window gives an introduction to sensors and how they respond to change. Up to four section displays can be open at any one time and they can be any combination of sensors and displays.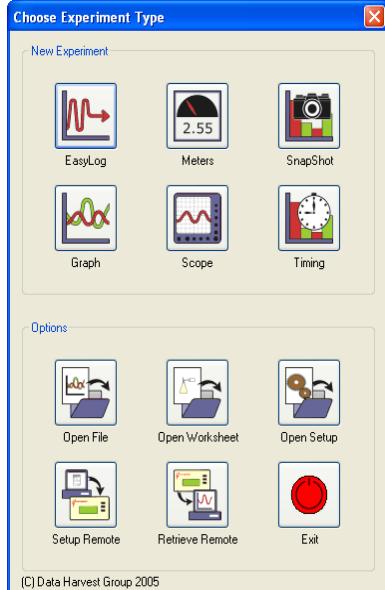

The Home page at Level 2 and 3

**Note:** If the interface/data logger detected cannot use an option it will appear grayed out e.g. the Setup and Retrieve Remote options will only be available for data loggers which support remote data collection.

## English

The program will open with live data being displayed from a sensor in a Numeric window. Click on the appropriate picture icon on the Menu bar to open other section displays i.e. Numeric, Gauge, Color Change or Bar. To alter the sensor displayed in a section, **right** click in the individual section window and select the sensor from the list.

Click on the Start/Stop icon  to stop recording.

### Snapshot

The Graph window will open pre-set to record sensor values on demand (**manual sampling**) with a bar graph display.

<input checked="" type="checkbox"/>
<input checked="" type="checkbox"/>
<input checked="" type="checkbox"/>
Level 1, 2, 3

Click on the Start/Stop icon  to start recording. Click in the graph area to record the current value for each active sensor. Click on the Start/Stop icon  to stop data being recorded.

### Graph

Graph is used to record sensor values against a **chosen time span** as a line graph display. The window opens with a real time logging wizard from which the time span, interval between samples, start condition and triggers for the recording can be chosen. (The options available will depend on the level you have selected to work at). When your selection has been made, click on Finish.

<input checked="" type="checkbox"/>
<input checked="" type="checkbox"/>
<input checked="" type="checkbox"/>
Level 1, 2, 3

Click on the Start/Stop icon  . Recording will start when the start condition you have selected is met e.g. if None is the selected trigger then the recording will start as soon as the Start/Stop icon is clicked. The recording will stop when the selected time span has passed.

**Note:** If you want to stop a recording before it has run the full term, click on the Start/Stop icon. 

### Scope

The Graph window will open with a simple oscilloscope type display, ideal for demonstrating fast events where there are rapid changes e.g. ECG traces, sound waveform. Click on the Start/Stop icon. 

<input checked="" type="checkbox"/>
<input checked="" type="checkbox"/>
<input checked="" type="checkbox"/>
Level 2, 3

The data from the sensor will be redrawn continuously on the screen.


## Timing

Timing is used to allow students to study Time, Velocity, Acceleration, Momentum and Kinetic Energy relationships using switch-type (digital) sensors such as Light Gates. A recording wizard will open with these options available to select before recording data. Sensors are connected to inputs A & B. (The options available will depend on the level you have selected to work at). When the selection has been made, click on Finish.

✓
✓
✓
Level 1, 2, 3

Click on the **Start/Stop**  icon and a reading will be displayed when a change in signal is detected from the digital sensors.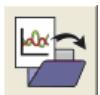

## Open File

Used to load a previously saved .ssl or .sid file. Locate the correct file and click on open.


✓
✓
✓
Level 1, 2, 3

## Open Worksheet

To select and open a worksheet and its accompanying experimental set up file.

✓
✓
✓
✓
Level 1, 2, 3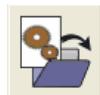

## Open Setup

To select and open an experimental set up file. This will preset the software for a specific operation.

✓
✓
x
Level 2, 3

## Set up Remote

Used to set up the available options for capturing data while disconnected from the computer (remote). These options can include the duration of a recording, the time between each sample and triggers. (The options available will depend on the level you have selected to work at).

✓
✓
x
Level 2, 3

## Retrieve Remote

Use this option to retrieve a data set stored in your data logger's memory.

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